



# INSTRUCTION MANUAL

**505857**

Biological Microscope

[www.wpiinc.com](http://www.wpiinc.com)

082724

---

CONTENTS

ABOUT THIS MANUAL ..... 1

INTRODUCTION ..... 1

    Parts List..... 1

    Unpacking..... 2

    Notes and Warnings..... 3

INSTRUMENT DESCRIPTION ..... 4

    Setup ..... 6

TROUBLESHOOTING ..... 13

WARRANTY ..... 18

    Claims and Returns ..... 18

    Repairs..... 18

*Copyright © 2024 by World Precision Instruments. All rights reserved. No part of this publication may be reproduced or translated into any language, in any form, without prior written permission of World Precision Instruments, Inc.*



## ABOUT THIS MANUAL

The following symbols are used in this guide:



This symbol indicates a CAUTION. Cautions warn against actions that can cause damage to equipment. Please read these carefully.



This symbol indicates a WARNING. Warnings alert you to actions that can cause personal injury or pose a physical threat. Please read these carefully.

NOTES and TIPS contain helpful information

## INTRODUCTION

### Parts List

After unpacking, verify that there is no visible damage to the sensor. Verify that all items are included:

(4) Plan Achromat Objective Lenses: 4x, 10x, 40xR, 100xR (oil immersion)\*

(1) Pair of Eyepieces, Widefield High Eyepoint Focusing 10x eyepieces/FN 20mm

(1) Trinocular Siedentopf Viewing Head, 30° inclined binocular eyepiece with interpupillary adjustment 47-75mm

(1) Vertical Trinocular Tube for camera integration

(1) Specimen Stage, 150mm x 140mm mechanical stage

(1) Illumination, Fixed Koehler illuminator, 3W LED with variable rheostat control and on/off switch, 110~240V power supply

(1) Camera, HD 1080p camera, live resolution 1920x1080 pixels, 2MP image capture to USB flash drive (drive not included), iLive image 1080p/60 fps (HDMI connection to 13" HD monitor)

(1) 13" LCD HDMI Monitor, 1080p with 110~240v power cord and HDMI cable

(1) Microscope Base with Cord and Accessory Storage Compartment (located on the back of the microscope frame)

(1) Instruction Manual is available online at [www.wpiinc.com/manuals](http://www.wpiinc.com/manuals).

\*All microscope objectives are parfocussed, parcentered and color-coded

---

## Unpacking

Upon receipt of this instrument, make a thorough inspection of the contents and check for possible damage. Missing cartons or obvious damage to cartons should be noted on the delivery receipt before signing. Concealed damage should be reported at once to the carrier and an inspection requested. Please read the section entitled "Claims and Returns" on page 18 of this manual. Please contact WPI Customer Service if any parts are missing at 941.371.1003 or [customerservice@wpiinc.com](mailto:customerservice@wpiinc.com).

**Returns:** Do not return any goods to WPI without obtaining prior approval (RMA # required) and instructions from WPI's Returns Department. Goods returned (unauthorized) by collect freight may be refused. If a return shipment is necessary, use the original container, if possible. If the original container is not available, use a suitable substitute that is rigid and of adequate size. Wrap the instrument in paper or plastic surrounded with at least 100mm (four inches) of shock absorbing material. For further details, please read the section entitled "Claims and Returns" on page 18 of this manual.

## Notes and Warnings



Do not shake or drop the microscope.




Do not expose the microscope to direct sun, high temperatures, high humidity, dust, or damp environments. Use a flat work surface.



When moving the microscope use both hands, holding the handle at the back (1) and the frame at the front (2) as shown in the image at right.




Set the power switch to off "O" before replacing a bulb or fuse, and wait until the lamp is cool. The power switch is located on the side of the microscope.

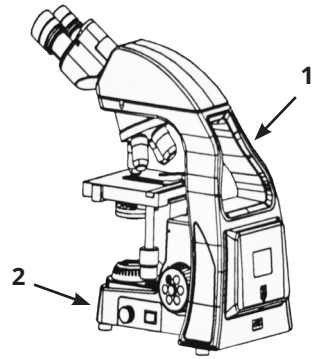


Voltage range of 100~240V (47-63Hz) is supported.




Turn the rheostat control knob (1) to the minimum and turn the power switch to the on "I" position. Rotate the rheostat knob to increase or decrease the brightness of the illumination.

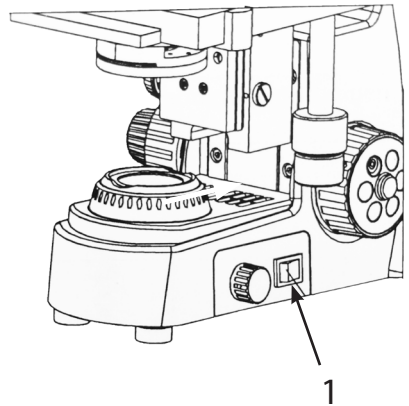

# INSTRUMENT DESCRIPTION

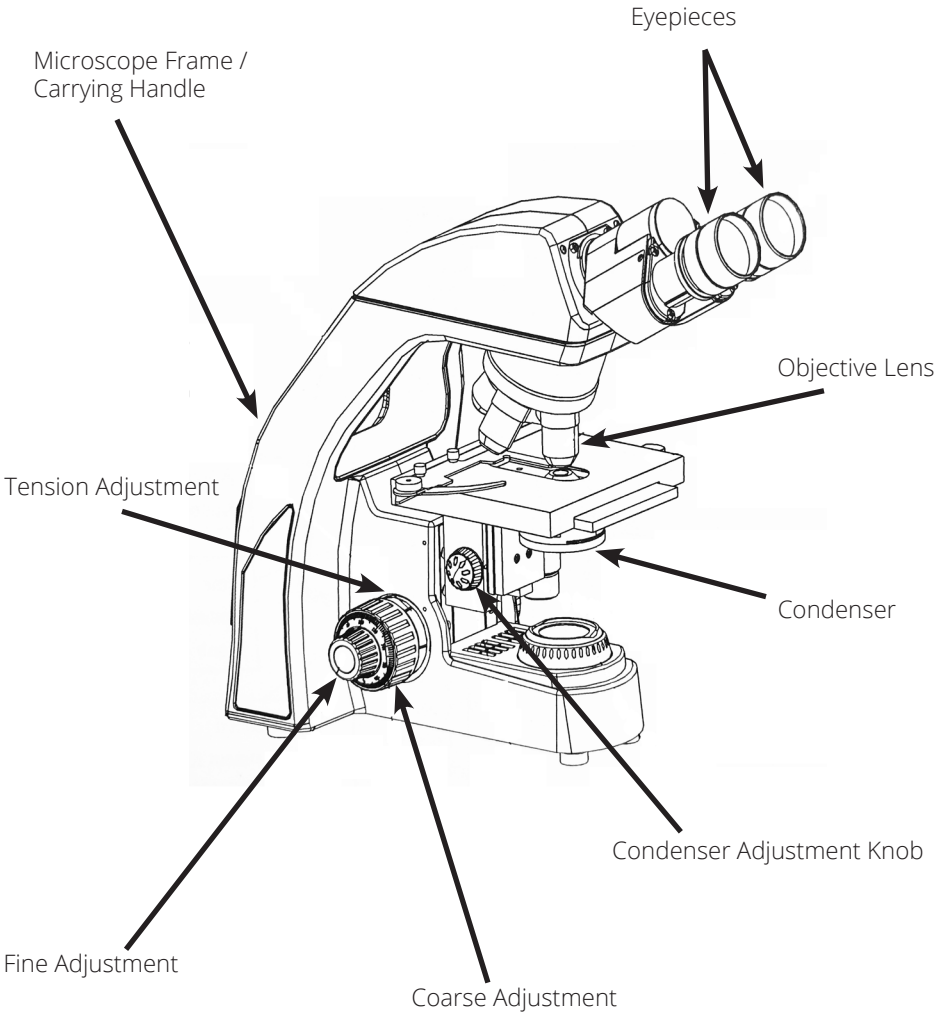




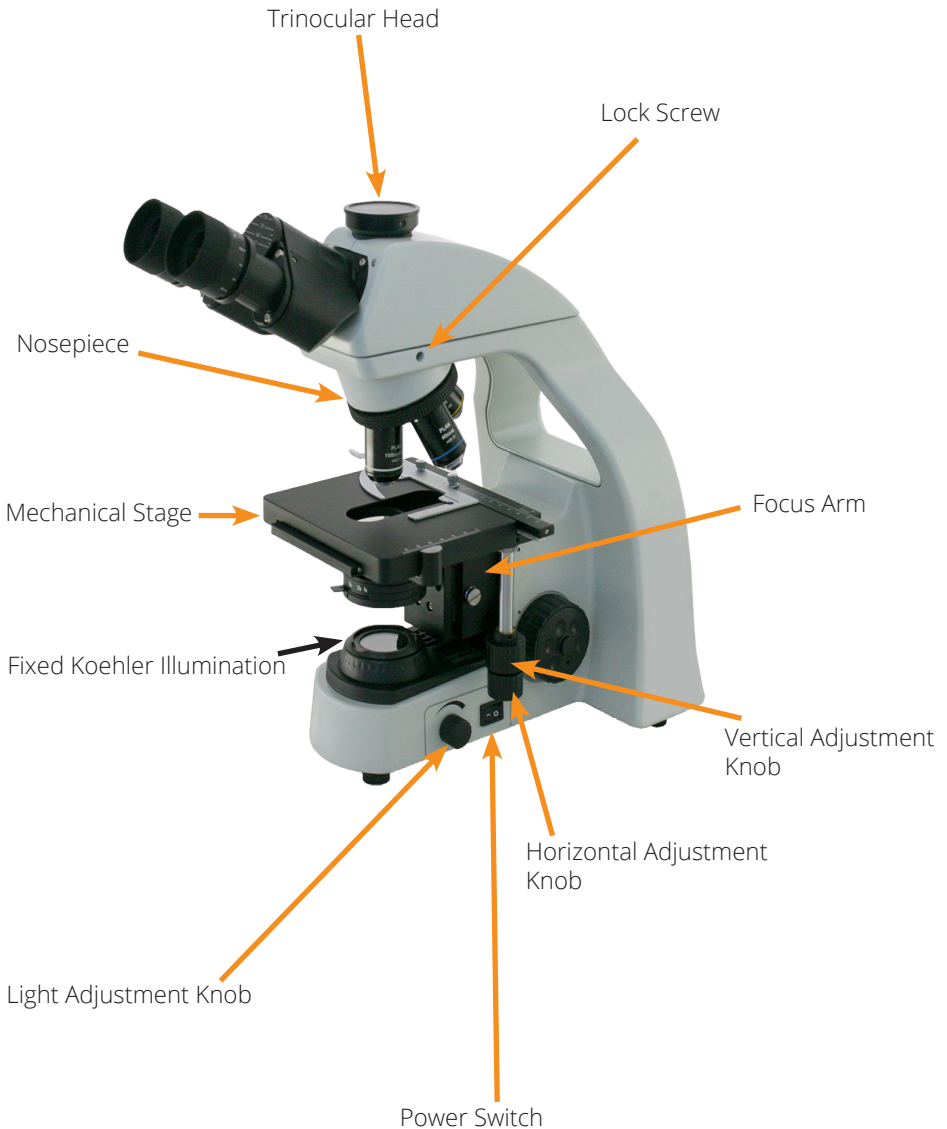

# Setup

Follow the numbers below when assembling the microscope.

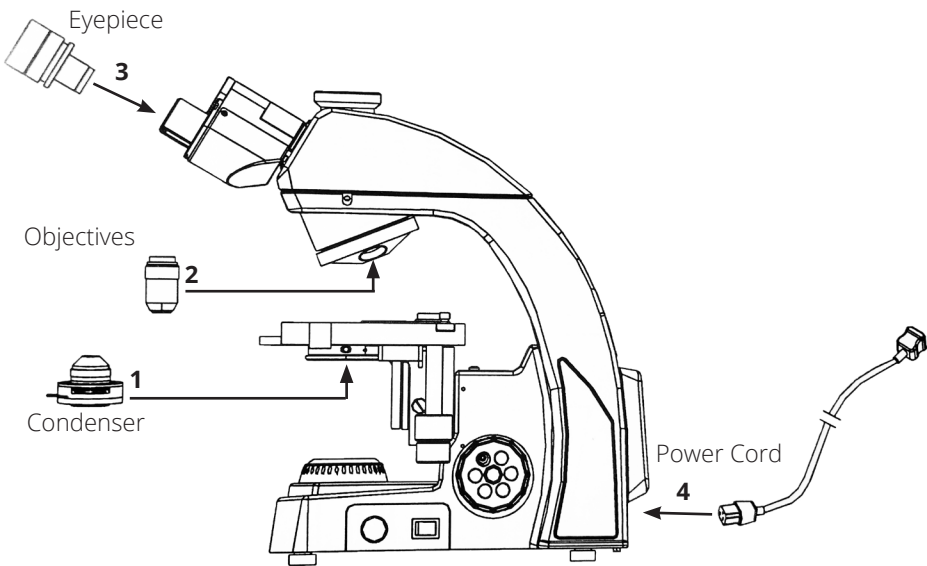




## Assemble the Condenser:

Rotate the coarse focus knob (1) to raise the stage to its highest position.


Fully loosen the condenser lock-screw (2).

Insert the condenser into the hole of base according to the arrowhead pointed (3), until the condenser is aligned with the base, and rotate the condenser to make the handle face forward. Tighten the condenser lock-screw (2) of condenser, then raise the condenser with the up-down knob to the highest position.

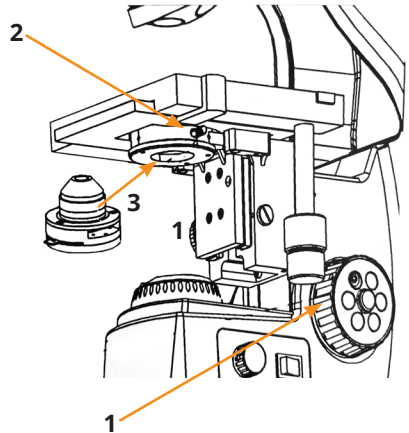

## Assemble the Objective:

Rotate the coarse focusing knob (1) to lower the stage to the lowest position. Then install the objectives into the nosepiece (2) from the lowest magnification to the highest in a clockwise direction.

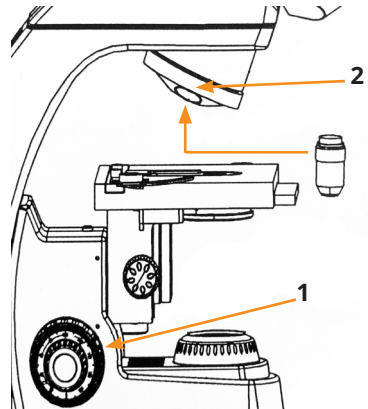




### Assemble the Eyepiece:

Take off the eyepiece tube cover (1).

Insert the eyepiece (2) into the eyepiece tube, until the eyepiece is flush with the eyepiece tube.

Tighten the eyepiece with M2.5 inner hexagon lock-screw (3).

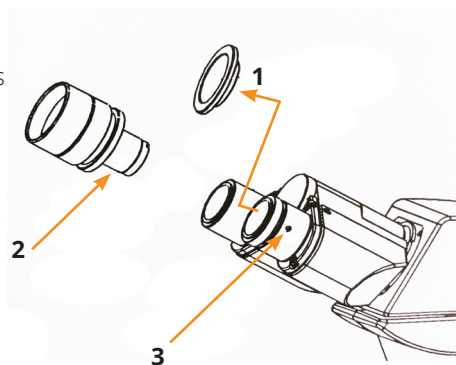


### Assemble the Eyepiece:

Take off the eyepiece tube cover (1).

Insert the eyepiece (2) into the eyepiece tube, until the eyepiece is flush with the eyepiece tube.


Tighten the eyepiece with M2.5 inner hexagon lock-screw (3).

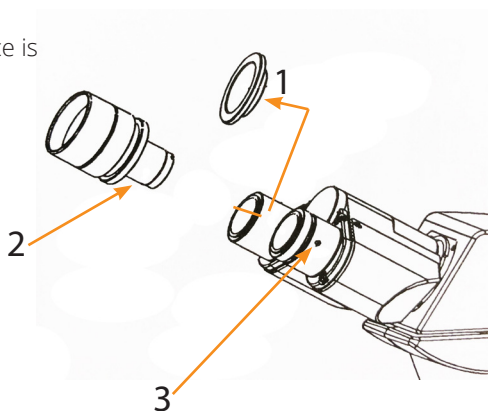




## Turning the Lamp ON:

Set the power switch to the ON position “-” (1). Adjust the illumination knob rotating clockwise to increase the light intensity and rotating counterclockwise to decrease the light intensity.

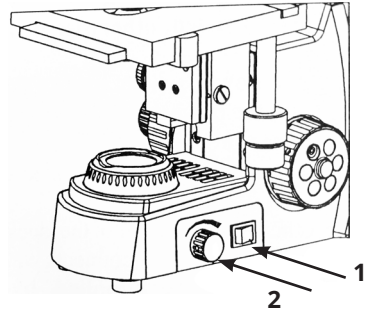


## Placing the Specimen Slide:

Push the spring lever (1) of the specimen holder backwards. Place the specimen slide (2) into the clip, then slowly move the specimen holder lever back into place. This will secure the specimen slide. Rotate the X and Y axis knobs (3) on the stage until the center of the specimen slide is aligned with the objective.

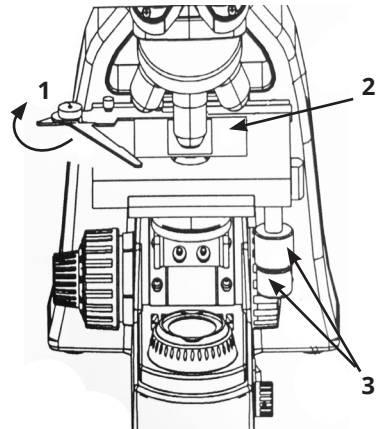

## Adjust the Focus:


Rotate the condenser adjustment knob (1) raising it to the highest position. Move the 4x objective into the optical path. Looking through the eyepieces, rotate the coarse focusing knob (2) until the specimen appears in the field of view and begins to come into focus. Rotate the fine focusing knob (3) to bring the specimen into crisp focus.

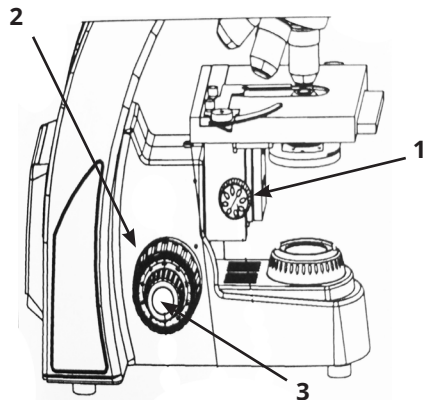




### Adjust the Focusing Tension:

If it becomes hard to move the coarse focus knob, or the stage drifts and the sample falls out of focus the tension needs to be adjusted. Tighten the tension by adjusting the tension adjustment ring (1) in the direction of the arrow shown at right. Loosen the tension by rotating the tension adjustment ring in the opposite direction.

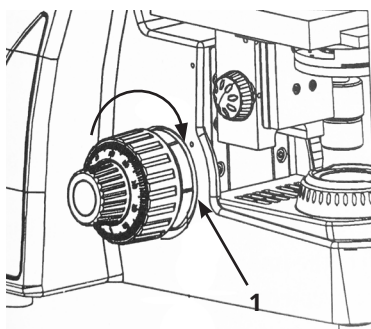


### Adjusting the Interpupillary Distance:

When observing with both eyes, adjust the eyepieces (2) to fit your personal interpupillary distance. The scale between the eyepieces (1) notes the interpupillary distance setting. It's a good idea to note this number when working in a setting where multiple people are using the same instrument. This allows you to easily reset your interpupillary distance.

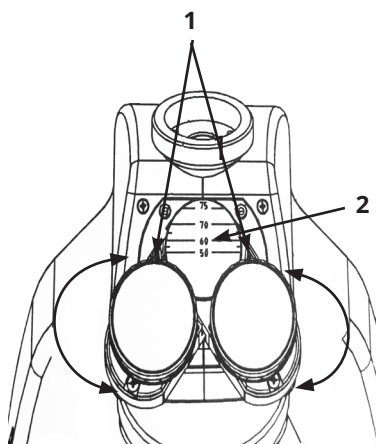

### Aperture Diaphragm:


The aperture diaphragm dictates the numerical aperture (NA) of the illumination. The best resolution, contrast, and depth of field is obtained when the NA of the illumination matches with the NA of the objective. Rotate the diaphragm adjustment ring (1) so the pointer matches up with the inscription of the objective being used.

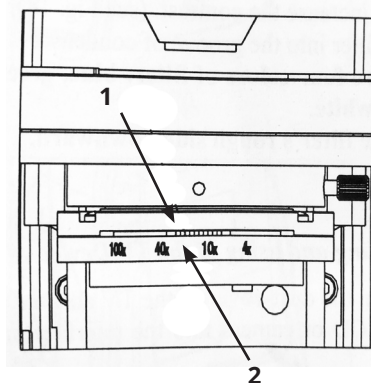




## Using the 100x Immersion Oil Objective:

Start with the 4x objective and focus on the sample. Place a drop of immersion oil (1) on the specimen. Rotate the nosepiece counterclockwise so the 100x objective is in the light path and the oil seals the space between the objective and the cover slip. Use the fine focus knob to obtain a clear image. Make sure there are no air bubbles in the oil. If there are, rotate the nosepiece slightly to remove them. After use, wipe the lens with a tissue moistened with microscope cleaning solution.

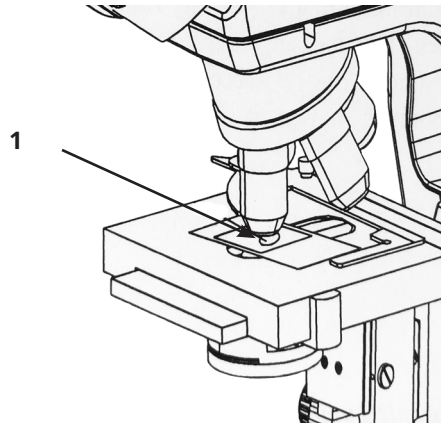


## Using Filters:

Filters can increase the contrast in microscopy images. The filter sits on top of the light as shown at the right. The 505857 accepts a 45mm diameter filter.

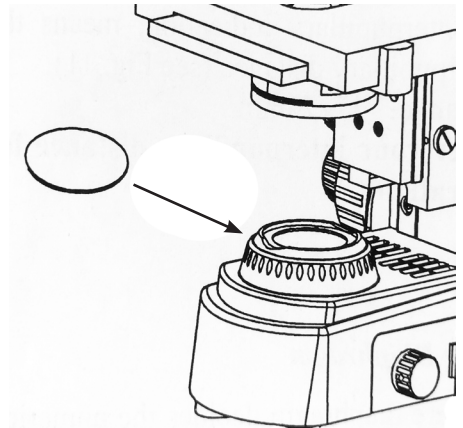

## C-Mount assembly:



Loosen the trinocular head set screw (1) and remove the trinocular port dust cap (2). Remove any dust cover caps from the c-mount (3). Insert the c-mount into the trinocular port as image shows at right and tighten the set screw (1). Connect the camera to the c-mount threads (4) and re-tighten the set screw. When the eyepieces are in focus, adjust the focus on the c-mount to parfocal the camera with the eyepieces.

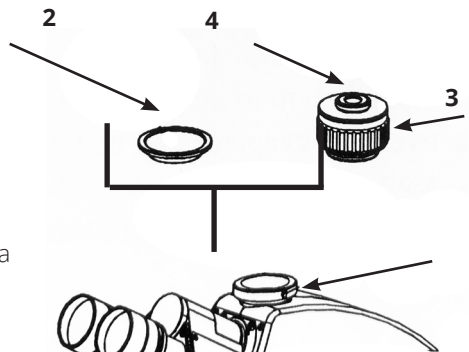



### Using the Storage Box:

When the microscope is not in use, the power cord can be placed in the storage box. Press the position (1) of the storage box cover with thumb. When you hear a click, push up and put the power cord into the storage box. Close the back cover by pressing down on the cover.

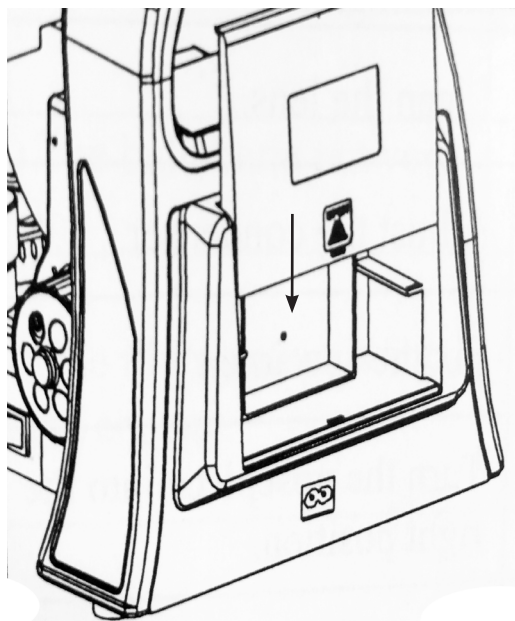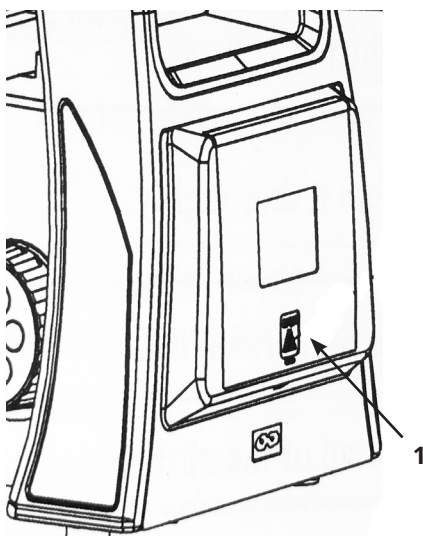



# TROUBLESHOOTING


TROUBLE	POSSIBLE CAUSE	SOLUTION
Field of view is dark even when light is on	The field diaphragm is not large enough.	Enlarge the field diaphragm
	Condenser is too low.	Adjust the position of condenser.
The edge of the field of view is dark or not even.	The nosepiece is not in the right position.	Turn the nosepiece into the right position.
	Stain or dust has accumulated on the lens such as condenser, objective, or eyepiece.	Clean the lens.
Stain or dust is observed in the field of view.	Stains have accumulated on the specimen.	Clean the specimen.
	Stains have accumulated on the lens.	Clean the lens.
Glare or hot spot on observed image.	Condenser is set too low.	Raise condenser.
	Aperture iris diaphragm is closed too much.	Open aperture iris diaphragm.
Unclear image	There's no cover glass on the specimen.	Add the cover glass.
	The cover glass is not standard.	Use a standard cover glass with thickness of 0.17mm.
	The slide is upside down.	Adjust the slide.
	The immersion oil has accumulated on the dry objective.	Clean the objective thoroughly.
	The immersion oil is not being used for oil objective.	Use immersion oil.
	Air bubble is in the immersion.	Get rid of the air bubbles.
	The aperture diaphragm is not opened correctly.	Adjust the aperture diaphragm.
	Stain or dust has accumulated on the eyepiece lens.	Clean the lens.
	The condenser is too low.	Adjust the condenser.

TROUBLE	POSSIBLE CAUSE	SOLUTIONN
One side of the field of view is dark or the image moves while in focus.	The specimen is not fixed.	Fix the specimen.
	The nosepiece is not in the right position.	Turn the nosepiece into the right position.
	Condenser centered incorrectly.	Center the condenser.
The eyes feel tired easily. The right field of view doesn't superpose with the left.	Interpupillary distance is wrong.	Adjust the interpupillary distance.
	Eyepieces for the left eye and right eye are different.	Use the same eyepiece.
Coarse focusing knob is too tight.	Tension knob is too loose.	Tighten it a little.
Stage declines itself, cannot stay on the focal plane when observing.	Tension knob is too loose.	Tighten it a little.
Coarse focusing knob can't rise.	The limit stop knob is locked.	Loosen the knob.
Coarse focusing knob can't decline.	The base of the condenser is too low.	Raise the base.
Cannot move the slide smoothly.	The slide is not fixed correctly.	Adjust it
	The movable specimen holder is not fixed properly.	Adjust it
The image moves obviously when touching the stage.	The stage is fastened incorrectly.	Fasten the stage correctly.
The LED light does not work.	No power supply.	Check the connection of the power cable.
	The LED bulb is not installed correctly.	Install it correctly.
	The LED bulb is burnt out.	Replace it.
The bulb burns out frequently.	The wrong bulb is being used.	Replace it with the correct bulb.
The illumination is not bright enough.	The wrong bulb is being used.	Replace it with the correct bulb.
	The illumination adjustment is not set in a bright position.	Adjust the illumination knob to intensify the light.



Before performing any of these steps the microscope **MUST** be turned off and unplugged.


With the microscope powered off and unplugged, gently lay the microscope down on a sturdy surface ensuring the carrying handle side is down.

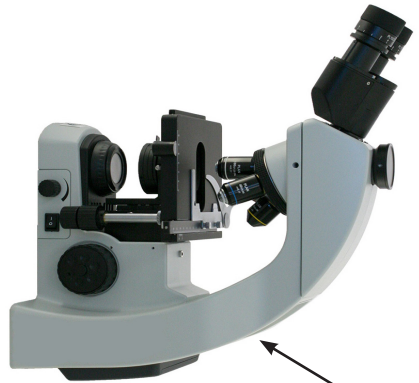


Gently lay microscope on carrying handle.



Locate the 4 rubber feet on the underside of the microscope. Remove the rubber feet exposing the 4 screws securing the bottom plate of the microscope.

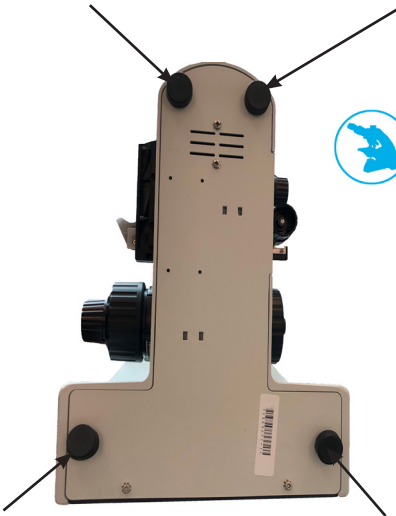


Using an Allen wrench remove the 4 screws securing the bottom plate of the microscope. With all 4 screws removed, gently remove the bottom plate exposing the internal electrical components of the microscope.

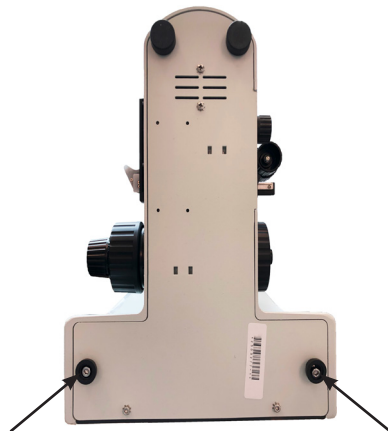



With all 4 screws removed, carefully remove the bottom plate exposing the internal components of the microscope.

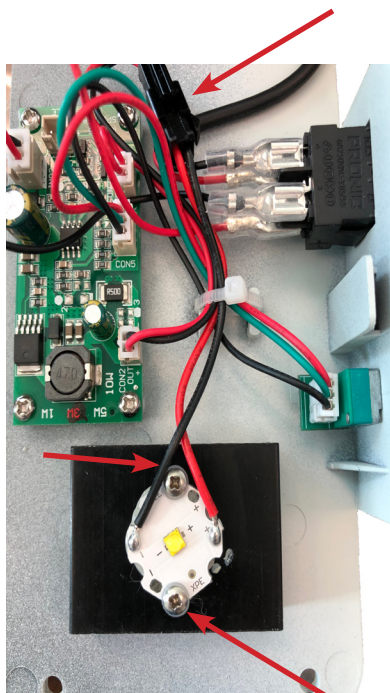

The electrical components are mounted onto the internal side of the bottom plate.



Locate the 2 screws securing the LED bulb in place. Using a screw driver, remove the 2 screws.



Locate the black wire connection. Using your thumb and forefinger, pinch the sides of the black wire plug connection and gently pull it from the wire housing.



Place the new LED bulb on to the LED bulb platform and attach using the small set screws that were previously removed.



LED replacement bulb part




Insert the black plastic wire plug back into the wire housing.



With the LED bulb replaced, position the bottom of the microscope plate back onto the microscope frame. Using a screw driver, secure the bottom plate with the 4 set screws. Place the 4 rubber feet over the top of the 4 set screws.



Return the microscope to its upright position. Plug the microscope back in, power on, and adjust the rheostat control as needed.

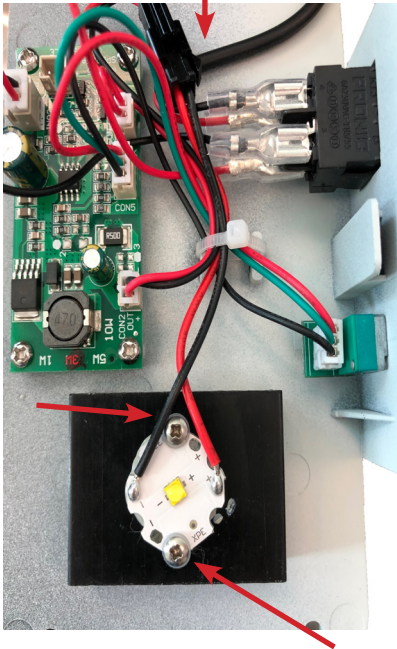

---

## WARRANTY

WPI (World Precision Instruments) warrants to the original purchaser that this equipment, including its components and parts, shall be free from defects in material and workmanship for a period of one year\* from the date of receipt. WPI's obligation under this warranty shall be limited to repair or replacement, at WPI's option, of the equipment or defective components or parts upon receipt thereof f.o.b. WPI, Sarasota, Florida U.S.A. Return of a repaired instrument shall be f.o.b. Sarasota.

The above warranty is contingent upon normal usage and does not cover products which have been modified without WPI's approval or which have been subjected to unusual physical or electrical stress or on which the original identification marks have been removed or altered. The above warranty will not apply if adjustment, repair or parts replacement is required because of accident, neglect, misuse, failure of electric power, air conditioning, humidity control, or causes other than normal and ordinary usage.

To the extent that any of its equipment is furnished by a manufacturer other than WPI, the foregoing warranty shall be applicable only to the extent of the warranty furnished by such other manufacturer. This warranty will not apply to appearance terms, such as knobs, handles, dials or the like.

WPI makes no warranty of any kind, express or implied or statutory, including without limitation any warranties of merchantability and/or fitness for a particular purpose. WPI shall not be liable for any damages, whether direct, indirect, special or consequential arising from a failure of this product to operate in the manner desired by the user. WPI shall not be liable for any damage to data or property that may be caused directly or indirectly by use of this product.

## Claims and Returns

Inspect all shipments upon receipt. Missing cartons or obvious damage to cartons should be noted on the delivery receipt before signing. Concealed loss or damage should be reported at once to the carrier and an inspection requested. All claims for shortage or damage must be made within ten (10) days after receipt of shipment. Claims for lost shipments must be made within thirty (30) days of receipt of invoice or other notification of shipment. Please save damaged or pilfered cartons until claim is settled. In some instances, photographic documentation may be required. Some items are time-sensitive; WPI assumes no extended warranty or any liability for use beyond the date specified on the container

Do not return any goods to us without obtaining prior approval and instructions from our Returns Department. Goods returned (unauthorized) by collect freight may be refused. Goods accepted for restocking will be exchanged or credited to your WPI account. Goods returned which were ordered by customers in error are subject to a 25% restocking charge. Equipment which was built as a special order cannot be returned.

## Repairs

Contact our Customer Service Department for assistance in the repair of apparatus. Do not return goods until instructions have been received. Returned items must be securely packed to prevent further damage in transit. The Customer is responsible for paying shipping expenses, including adequate insurance on all items returned for repairs. Identification of the item(s) by model number, name, as well as complete description of the difficulties experienced should be written on the repair purchase order and on a tag attached to the item.

*\* Electrodes, batteries and other consumable parts are warranted for 30 days only from the date on which the customer receives these items.*





WORLD  
PRECISION  
INSTRUMENTS