



WORLD
PRECISION
INSTRUMENTS

INSTRUCTION MANUAL

CLEANRAY-1000

Tabletop Ultraviolet Sterilizer



www.wpiinc.com

051220



WPI's portable ultraviolet (UV) sterilizer is perfect for quick disinfection. UV sanitation has proven to be highly effective against pathogens. It is safe and easy to use and comes with a one-year warranty.

This unit is ideal for use in homes, offices, washrooms, child care centers, hotels, beauty and nail salons, barber shops, gyms, spa centers, hospitals and dental offices. It works for many small items like phones, face masks, watches, jewelry, glasses, keys, pacifiers, pet accessories, toothbrushes, towels, beauty & hairdressing equipment, kitchen equipment, surgical instruments, tattoo instruments and many other daily necessities.


Features

- Rapid sterilization with high power UV light and ozone
- No added reagents required for sterilization
- Large sterilization capacity
- Easy access to the sterilization chamber
- Safety lock turns off the UV light before the cabinet opens
- Safe for users because dark blue lid blocks ultraviolet light
- 9000 hours average bulb life

Notes and Warnings



CAUTION: The unit uses an ultraviolet light. Do not look directly at the unshielded bulb when it is in operation.

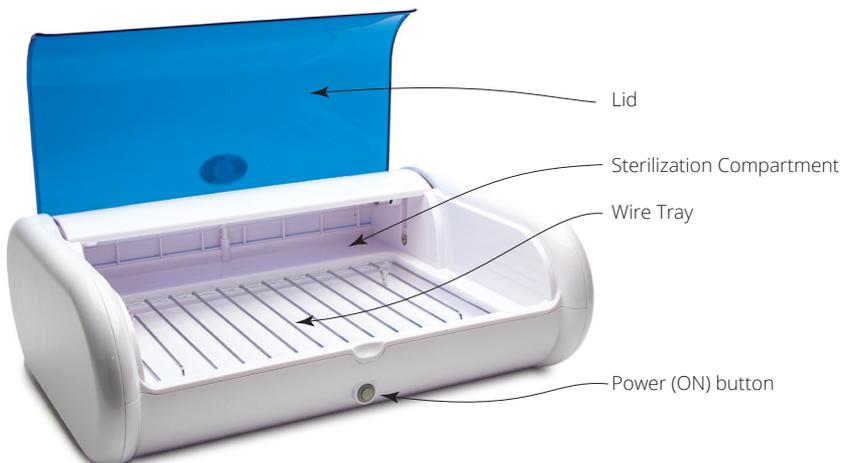

OPERATING INSTRUCTIONS

1. Plug in the unit to a wall AC outlet.
2. Arrange the items to be sterilized on the wire tray in a single layer.
3. Closed the lid.
4. Push the ON button located in the center of the front of the unit. A blue light will come on. Allow the light to run continuously for at least 15 minutes to sterilize the contents of the compartment. If you open the lid, the light automatically turns off.
5. Press the ON button again to turn the light off. Unplug the unit after use.

MAINTENANCE



CAUTION: Do not use acid, alkali, corrosive or toxic cleaners.

Using a pH neutral detergent solution, wipe the inside and outside with a damp cloth to remove dust and debris.

To change the UV lamp, open the cover, slide the bulb out and replace it with a new bulb. Dispose of the old bulb properly.

SPECIFICATIONS

This unit conforms to the following specifications:

Type	UV sterilizer
UVC Wavelength	254 nm
Safety	Interlock
Interior Volume	~ 6.5 L
Product Size	40 × 30 × 14 cm
Light Source	9 W, 3000 hour ave. bulb life
Voltage	110-120 V/220-240 V, 50/60 Hz
Plugs	EU, AU, US, UK
Material	ABS, AS cover, UV bulb
Weight	2 kg (4.4 lbs.)
Certification.....	CE

Warranty

WPI (World Precision Instruments) warrants to the original purchaser that this equipment, including its components and parts, shall be free from defects in material and workmanship for a period of 30 days from the date of receipt. WPI's obligation under this warranty shall be limited to repair or replacement, at WPI's option, of the equipment or defective components or parts upon receipt thereof f.o.b. WPI, Sarasota, Florida U.S.A. Return of a repaired instrument shall be f.o.b. Sarasota.

The above warranty is contingent upon normal usage and does not cover products which have been modified without WPI's approval and/or which have been subjected to unusual physical or electrical stress or on which the original identification marks have been removed or altered. The above warranty will not apply if adjustment, repair or parts replacement is required because of accident, neglect, misuse, failure of electric power, air conditioning, humidity control, or causes other than normal and ordinary usage.

To the extent that any of its equipment is furnished by a manufacturer other than WPI, the foregoing warranty shall be applicable only to the extent of the warranty furnished by such other manufacturer. This warranty will not apply to appearance terms, such as knobs, handles, dials or the like.

WPI makes no warranty of any kind, express or implied or statutory, including without limitation any warranties of merchantability and/or fitness for a particular purpose. WPI shall not be liable for any damages, whether direct, indirect, special or consequential arising from a failure of this product to operate in the manner desired by the user. WPI shall not be liable for any damage to data or property that may be caused directly or indirectly by use of this product.

Claims and Returns

- Inspect all shipments upon receipt. Missing cartons or obvious damage to cartons should be noted on the delivery receipt before signing. Concealed loss or damage should be reported at once to the carrier and an inspection requested. All claims for shortage or damage must be made within 10 days after receipt of shipment. Claims for lost shipments must be made within 30 days of invoice or other notification of shipment.

Please save damaged or pilfered cartons until claim settles. In some instances, photographic documentation may be required. Some items are time sensitive; WPI assumes no extended warranty or any liability for use beyond the date specified on the container.

- WPI cannot be held responsible for items damaged in shipment en route to us. Please enclose merchandise in its original shipping container to avoid damage from handling. We recommend that you insure merchandise when shipping. The customer is responsible for paying shipping expenses including adequate insurance on all items returned.
- Do not return any goods to WPI without obtaining prior approval and instructions (RMA#) from our returns department. Goods returned unauthorized or by collect freight may be refused. The RMA# must be clearly displayed on the outside of the box, or the package will not be accepted. Please contact the RMA department for a request form.
- Goods returned for repair must be reasonably clean and free of hazardous materials.
- A handling fee is charged for goods returned for exchange or credit. This fee may add up to 25% of the sale price depending on the condition of the item. Goods ordered in error are also subject to the handling fee.
- Equipment which was built as a special order cannot be returned.
- Always refer to the RMA# when contacting WPI to obtain a status of your returned item.
- For any other issues regarding a claim or return, please contact the RMA department

Warning: This equipment is not designed or intended for use on humans.

WORLD PRECISION INSTRUMENTS

USA: 175 Sarasota Center Boulevard, Sarasota FL 34240-9258 USA

Tel: (941) 371-1003 • **Fax:** (941) 377-5428 • wpi@wpiinc.com • www.wpiinc.com

UK: 1 Hunting Gate, Hitchin, Hertfordshire SG4 0TJ England • 44 (0)1462 424700 • wpiuk@wpi-europe.com

Germany: Pfingstweide 16, D-61169 Friedberg (Hessen) • Tel: +49 (0)6031 67708-0 • Fax: +49 (0)6031 67708-80 • wvide@wpi-europe.com

China & Hong Kong: Rm 27A, No 8 Donfang Rd., Pudong New District, Shanghai PRC • +86 68885517 • ChinaSales@china.wpiinc.com

Brazil: Conselheiro Nabias, 756 sala2611, Santos-São Paulo 11045-002 Brazil • info@brazil.wpiinc.com