

Kima-Connect: Tubing set for Kima pump

Kima Connect tubing set for Kima pump contains:

- 2 pump priming cables
- 1 pump inlet cable
- 1 pump outlet cable
- 1 biochip outlet cable

To connect the tubing to the Kima pump, follow the steps illustrated below.



- 1.** Remove cover from pump. Connect pump priming cables to inlet and outlet of pump and prime with solution (e.g. deionised water or PBS) using the syringe.



- 2.** Connect inlet cable as shown.



- 3.** Replace cover on pump. Connect pump outlet cable as shown.



- 4.** Connect biochip outlet cable as shown.

