


## Mirus-Connect: Tubing set for Mirus Evo Nanopump

Mirus Connect tubing set for Mirus Evo Nanopump contains:

- 1 pump inlet cable
- 1 pump outlet cable
- 1 pump-to-MultiFlow8 cable

To connect the tubing to the Mirus Evo Nanopump, follow the steps illustrated below.

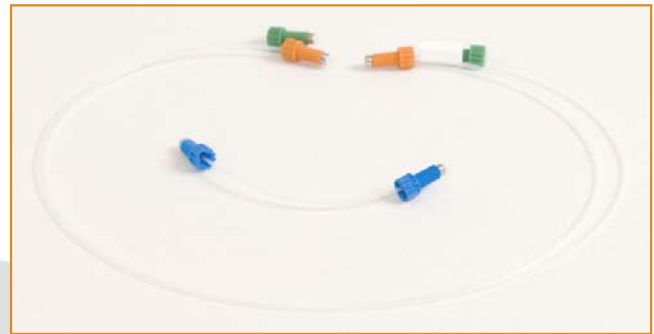

- 1.** Connect inlet cable from flask to pump as shown.



- 2.** Connect outlet cable as shown.



- 3.** Connect cable to MultiFlow8 as shown.

